

AMI 01 AMINES

Analysis of Amines on BPX5

KEYWORDS:


Amines
Aromatic Amines
Diamines

Notes: BPX5 is a chemically inert column ideal for aromatic amine analysis

Part No.: 054113
Phase: BPX5, 0.25 μ m
Column: 25m x 0.22mm I.D.
Initial Temp.: 40°C, 1 min
Rate: 8°C/min
Final Temp.: 360°C
Detector: FID, 380°C
Injector: Split
Carrier Gas: H₂, 12 psi

Components (5 ng)

1. Decylamine
2. 1, 9-Diaminononane
3. 4-Nitroaniline
4. 3-Nitroaniline
5. 2-Nitroaniline

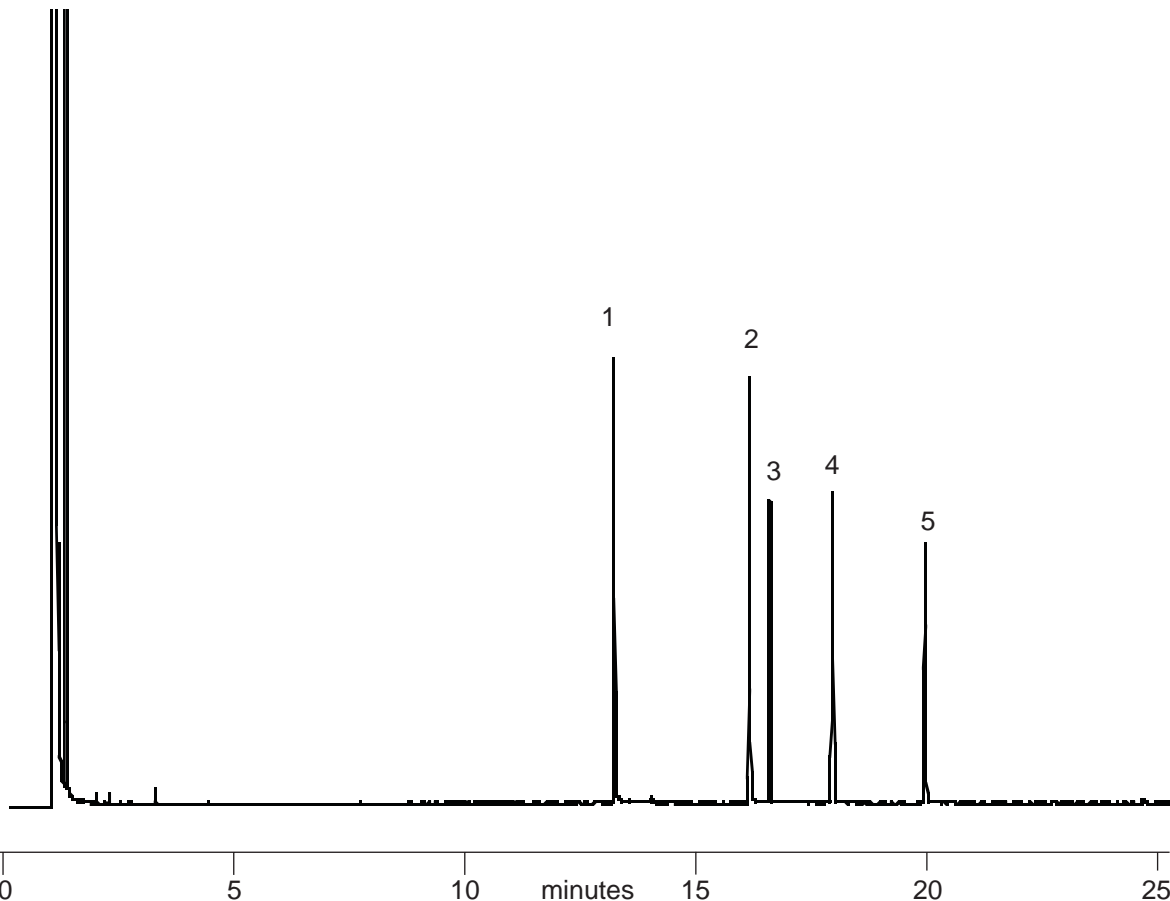

visit www.sge.com or contact our technical Customer Support Team on: techsupport@sge.com

SGE International Pty. Ltd.
email: support@sge.com

SGE Incorporated (USA)
email: usa@sge.com

SGE (France) Sarl
email: france@sge.com

SGE Deutschland GmbH
email: germany@sge.com

SGE Japan Inc.
email: sgejpn@itjit.ne.jp

SGE Australia Pty. Ltd.
email: sales@sge.com

SGE Europe Ltd. (UK)
email: uk@sge.com

SGE (Italia) Srl.
email: sge.italia@tin.it

SGE China Service Centre
email: sge@bj.col.com.cn

SGE India
email: sgeindia@bom4.vsnl.net.in