

FOO 23
FOOD

Free Fatty Acid Analysis on BP21

KEYWORDS:


Free Fatty Acids

Volatile Fatty Acids

Part No.: 054474
Phase: BP21, 0.5µm Film
Column: 25m x 0.53mm ID
Initial Temp.: 90°C, 1min
Rate 1: 8°C/min
Final Temp.: 160°C
Detector: FID, 300°C
Injector Mode: Split, 100:1
Carrier Gas: Helium, 5ml/min
Injection Volume: 1µL

Components

1. Acetic Acid
2. Propionic Acid
3. Isobutyric Acid
4. Butyric Acid
5. Isovaleric Acid
6. Valeric Acid
7. Unknown
8. Isocaproic Acid

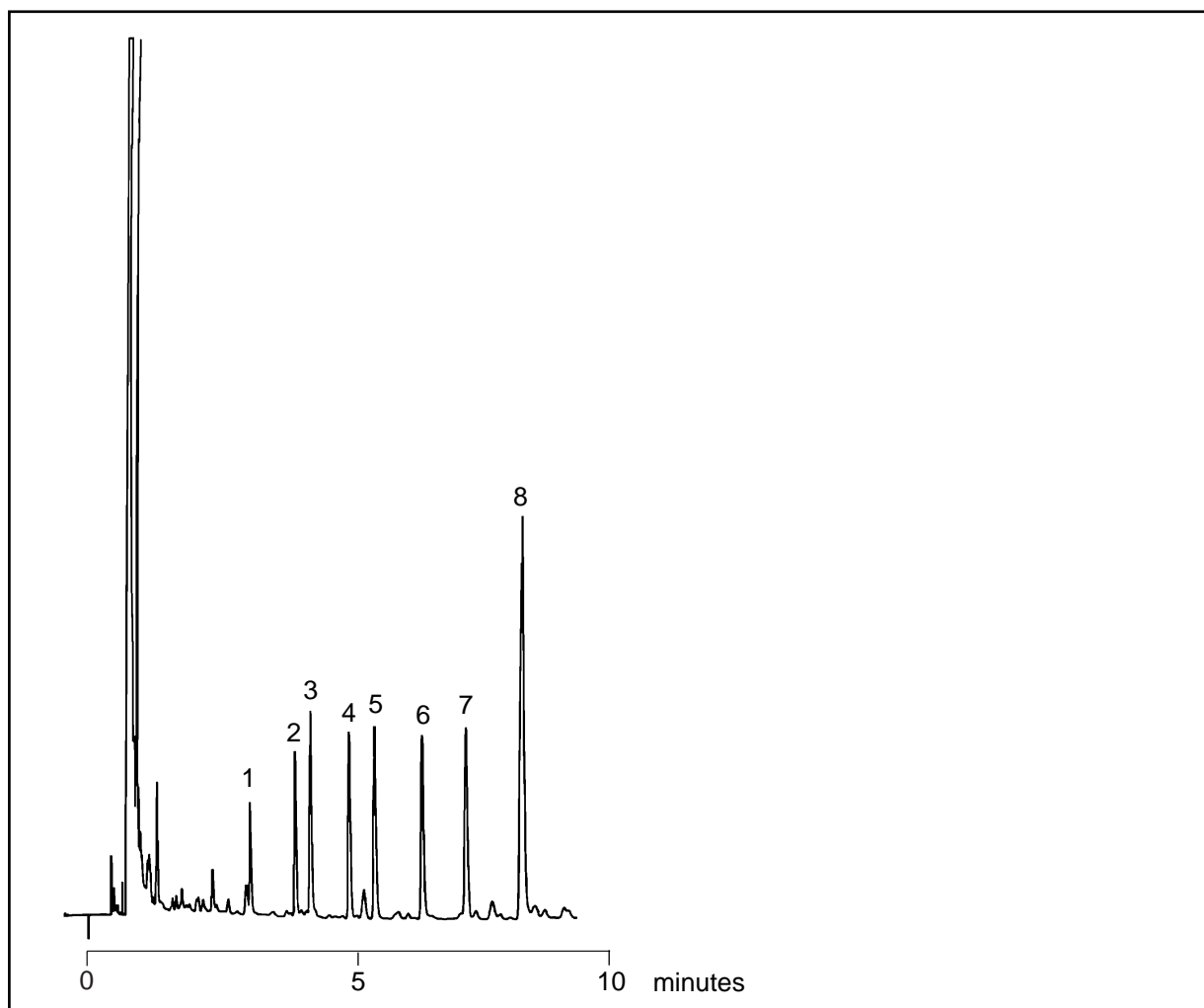

visit www.sge.com or contact our technical Customer Support Team on: techsupport@sge.com

SGE International Pty. Ltd.
email: support@sge.com

SGE Incorporated (USA)
email: usa@sge.com

SGE (France) Sarl
email: france@sge.com

SGE Deutschland GmbH
email: germany@sge.com

SGE Japan Inc.
email: sgejpn@itjit.ne.jp

SGE Australia Pty. Ltd.
email: sales@sge.com

SGE Europe Ltd. (UK)
email: uk@sge.com

SGE (Italia) Srl.
email: sge.italia@tin.it

SGE China Service Centre
email: sge@bj.col.com.cn

SGE India
email: sgeindia@bom4.vsnl.net.in