


GC COLUMN REPORT

Guide for Interpretation

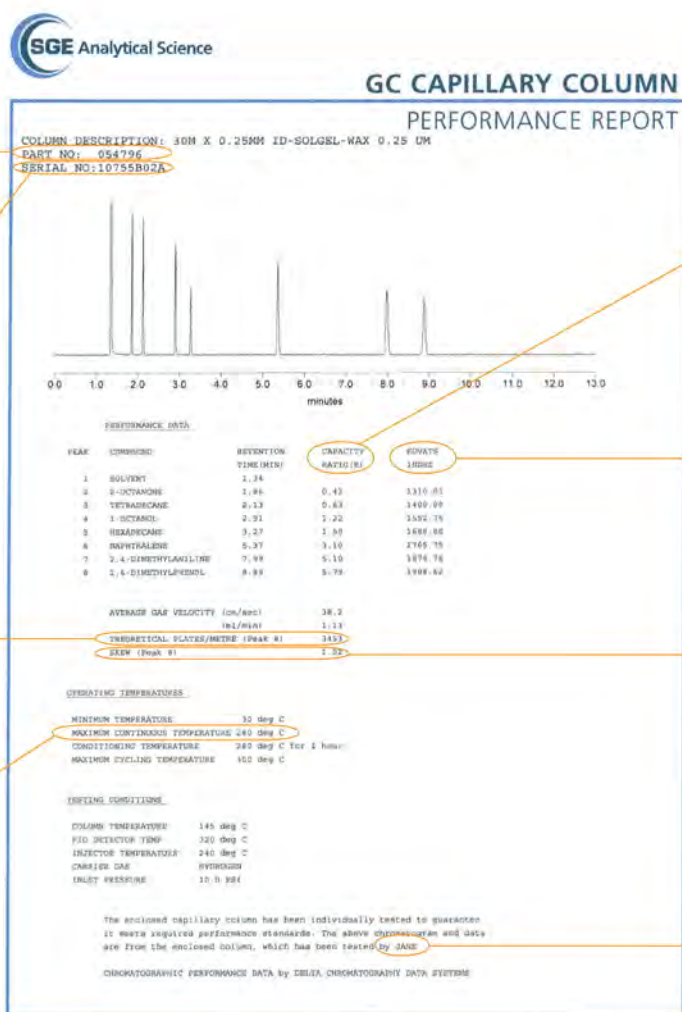

Part No.:
Re-order information.

Serial No. Column Traceability:
Every SGE GC column is traceable back to its manufacture.

Measure of Theoretical Plates/meter:
This is a measure of the efficiency of the column.

Maximum continuous temperature:
This is the maximum recommended temperature for the column. Higher temperatures can be used, but this will reduce column lifetime.

Capacity Ratio:
This is a measure of the film thickness.

Kovats Index:
Describes the retention behavior of a compound relative to that of straight chain hydrocarbons. Especially important for more polar columns.

Skew:
This is a measure of the degree of tailing (1.0 = Perfect).

Column Tester:
Every individual SGE column is tested to ensure quality.